

TOTAL PROTEIN STAIN FOR MEMBRANES

Amido Black

1. Stain filter for 1' in 0.1% amido black 10B in 25% isopropanol, 10% acetic acid.
2. Destain 2 x 15' in 25% isopropanol, 10% acetic acid.

Ponceau S (temporary stain that can be used prior to immunoreaction)

1. Stain filter for 1' in 0.2% Ponceau S in 3% TCA (trichloroacetic acid).
2. Wash filter with distilled H₂O until water runs clear and background staining is reduced.